
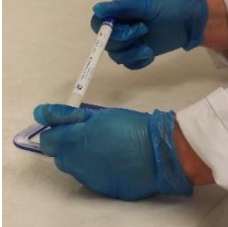
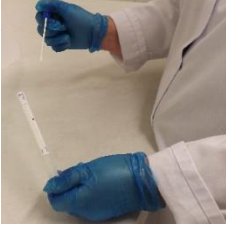
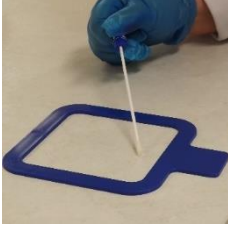

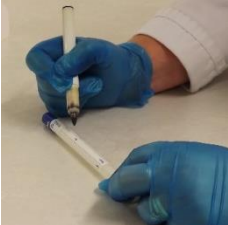


# INSTRUCTION FOR USE

## PreMoist Swab Kits

All samples should be kept cool until testing is performed.

<p><b>STEP 1</b></p> <p>Identify the site to be sampled.</p> 	<p><b>STEP 2</b></p> <p>Hold the swab with both hands and pull the swab cap away from the tube, breaking the security seal in the process.</p> 
<p><b>STEP 3</b></p> <p>Remove swab using the cap as a holder. Do not touch the swab tip or shaft and make sure that it does not touch any surface before sampling.</p> 	<p><b>STEP 4</b></p> <p>Keep the open transport tube with solution within reach, sample designated area using a template if possible. If the area permits, swab in 3 planes (horizontal, vertical and diagonal) whilst rotating.</p> 
<p><b>STEP 5</b></p> <p>When sampling is completed insert swab into the tube and push cap on tightly.</p> 	<p><b>STEP 6</b></p> <p>Write sample details on label, store in a cool (4°C -8°C), dry area and send to the laboratory within 24 hours for analysis.</p> 

Have you considered?

Sterile Sampling Template - TS/15-T



**Technical Service**  
Consultants Ltd

[www.tscswabs.co.uk](http://www.tscswabs.co.uk)  
Tel: +44 (0)1706 620600

IFU\_PreMoist\_Swab\_V1 –  
21/03/2014